

## Torch Tube EzyGlide Cloth - Instructions For Use

1. Disassemble your demountable torch.
2. Use protective gloves. The EzyGlide Cloth contains chemicals which may irritate the skin.
3. Apply an even coating of lubricant to the ground or 'grey' area of the torch tube.
4. Stand the tube upright and allow to dry for approximately 2 minutes before re-assembling the torch.
5. Replace the EzyGlide Cloth in the airtight container to prevent evaporation of the lubricant.

**Note:** The lubricant is designed to protect and lubricate in extreme conditions. Lubrication is only necessary if the torch tubes are difficult to insert into the torch body. For example: after acid cleaning of the torch tubes or extreme heating in the cleaning process.