




# VEESPRAY FITTING MAINTENANCE

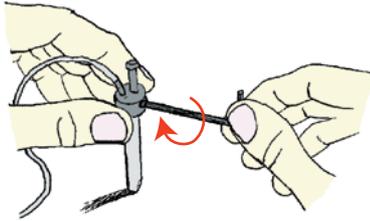


1. Holding the nebulizer in the left hand and leaning it vertically on a smooth flat surface, loosen the 2 grub screws with the right hand, using a small hex key.

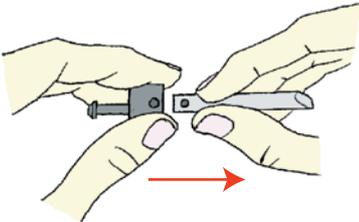

2. Take out the nebulizer body and the old sample tube.

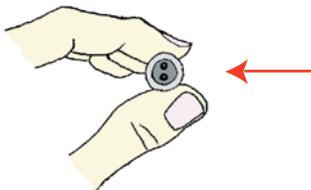


3. Check that the 2 o-rings in the nebulizer fitting are in place.

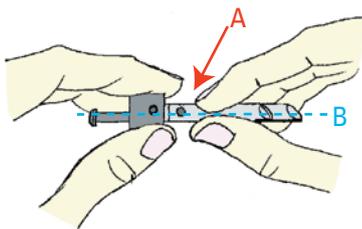


4 a. Replace the nebulizer body into the fitting, lining up the threaded holes with the indentations in the nebulizer body.

b. Line up the gas bore inside the nebulizer body with the gas arm.

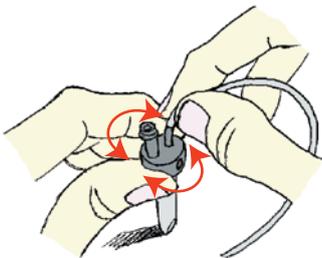

5. Holding the nebulizer in the left hand again, insert the new sample tube by twisting left to right and pushing it in gently, until it stops.

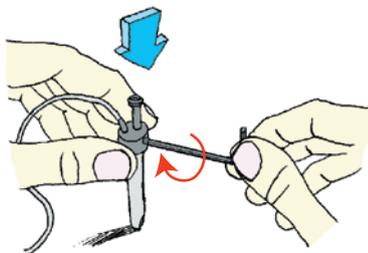

6. Ensure that the indentations in the nebulizer body line up with the threaded holes. Pushing down on the fitting, tighten the 2 grub screws firmly.

Note: Do not use excessive force when tightening.



**GLASS EXPANSION**

Quality By Design

**International**

Glass Expansion  
15 Batman Street  
West Melbourne VIC 3003  
Australia

Telephone: +61 3 9320 1111  
Toll Free (Aust): 1800 777 638  
Facsimile: +61 3 9320 1112  
Email: [enquiries@geicp.com](mailto:enquiries@geicp.com)

**Americas**

Glass Expansion  
4 Barlows Landing Road  
Unit 2A  
Pocasset, MA 02559, USA

Telephone: 508 563 1800  
Toll Free: 800 208 0097  
Facsimile: 508 563 1802  
Email: [geusa@geicp.com](mailto:geusa@geicp.com)

For your nearest Glass Expansion distributor, visit our website at [www.geicp.com](http://www.geicp.com)