


Instructions for Installing a Cyclonic Spray Chamber on the Varian Vista

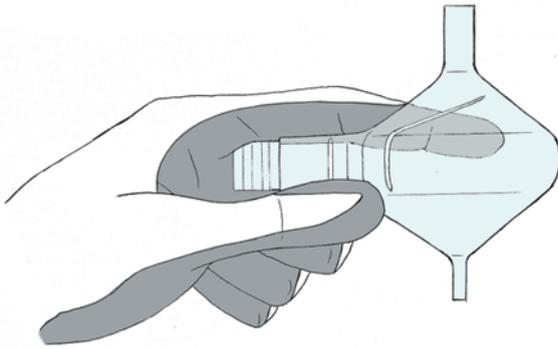

1. For your protection always wear safety gloves and safety glasses when handling glass.



2. Grip the spray chamber by the nebulizer arm (with the nebulizer removed).

3. Insert the aerosol outlet all the way into the upper slot on the mounting bracket.

4. Press down firmly to clip the drain outlet into the lower slot.

5. Connect the aerosol transfer tube and drain tube, and insert the nebulizer.

6. To remove the spray chamber, follow the above procedures in reverse.

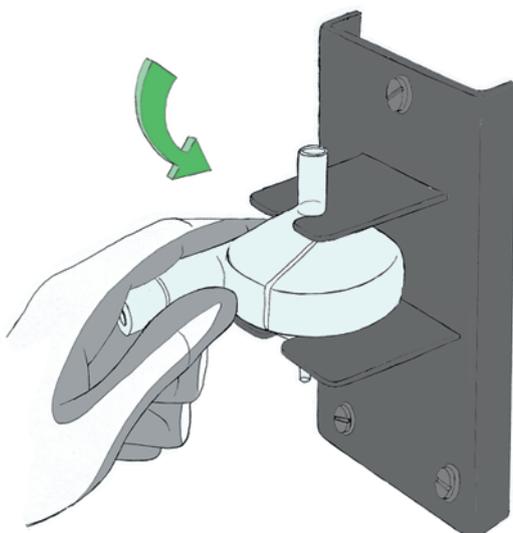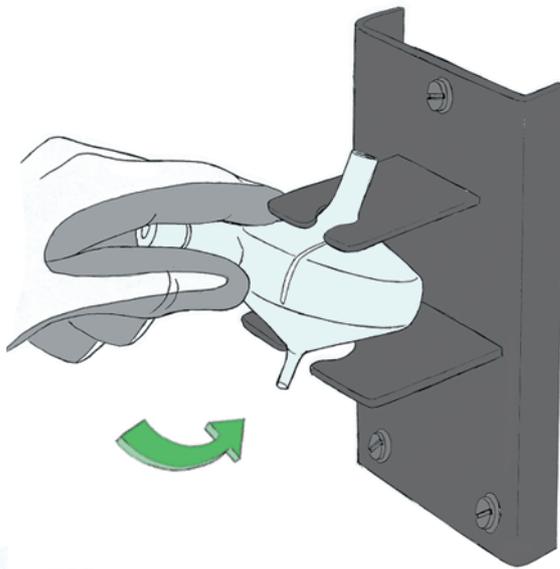

GLASS EXPANSION

Quality By Design

International

Glass Expansion
15 Batman Street
West Melbourne VIC 3003
Australia

Telephone: +61 3 9320 1111
Toll Free (Aust): 1800 777 638
Facsimile: +61 3 9320 1112
Email: enquiries@geicp.com

Americas

Glass Expansion
4 Barlows Landing Road
Unit 2A
Pocasset, MA 02559, USA

Telephone: 508 563 1800
Toll Free: 800 208 0097
Facsimile: 508 563 1802
Email: geusa@geicp.com

For your nearest Glass Expansion distributor, visit our website at www.geicp.com